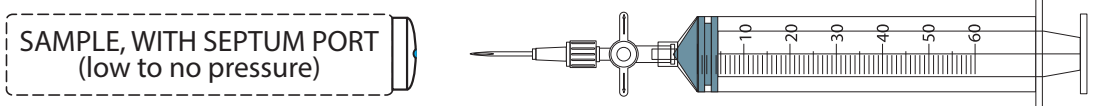
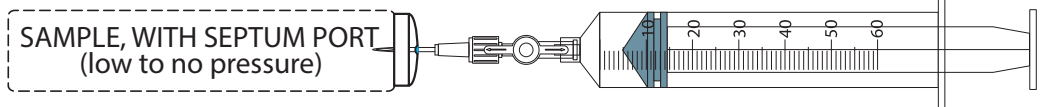
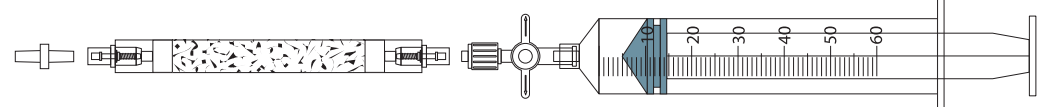
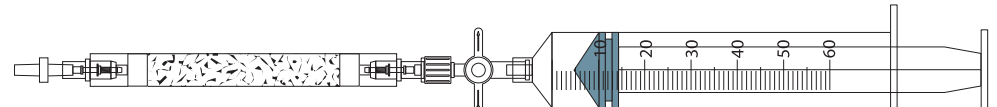
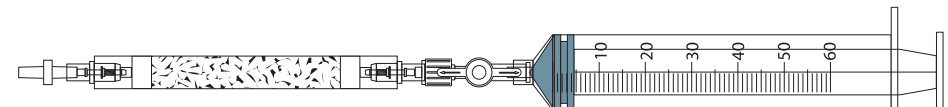
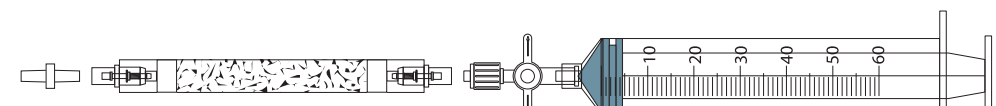
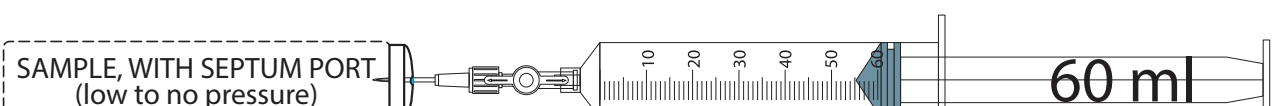
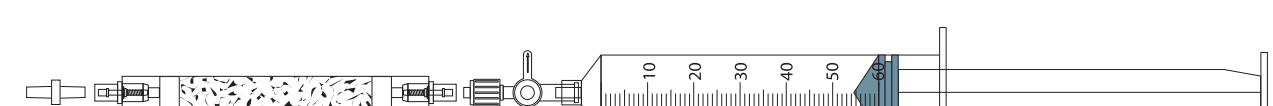
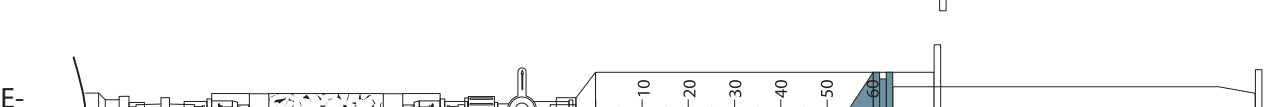


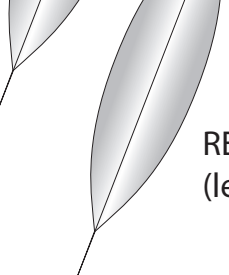


IsoScrubber™ Instructions

- 1.)  SAMPLE, WITH SEPTUM PORT
(low to no pressure)
- 2.) EXTRACT  SAMPLE, WITH SEPTUM PORT
(low to no pressure) **10 ml**
- 3.) DETACH NEEDLE 
- 4.) CONNECT SYRINGE & UNION 
- 5.) PURGE ISOSCRUBBER 
- 6.) DISCONNECT 
- 7.) EXTRACT  SAMPLE, WITH SEPTUM PORT
(low to no pressure) **60 ml**
- 8.) CLOSE VALVE-
DETACH NEEDLE 
- 9.) ATTACH SYRINGE-
ISOSCRUBBER-
UNION TO GAS BAG 
- 10.) FILL BAG 
- 11.) DETACH SYRINGE 
- 12.) REPEAT  REPEAT STEPS 7-10, TWO MORE TIMES
(leaving bag attached to IsoScrubber until 180 ml have been collected)